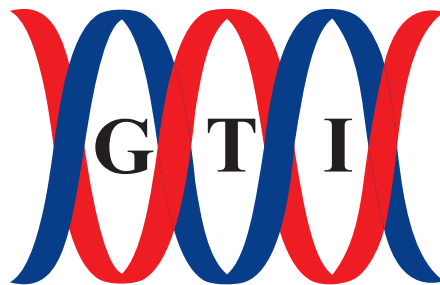

GTI-Microcentrifuge

Instruction Manual



Genetic Technology Instrumentation

Copyright © 2014



Operation

1. Turn on the unit.
2. Press the stop (red) button if the start light (green) is on.
3. Adjust the required time by turning the timer switch clockwise. Each click position represents 15 seconds. Time can be set from 15- 300 seconds. Setting at “~” position sets the timer to infinity.
4. Select speed (low/high).
5. Press the start (green) button to start the centrifuge.
6. Opening the lid automatically stops the centrifuge.
7. Centrifuge may be stopped prematurely by pressing the stop (red) button.
8. In case of power interruption the timer is reset to “off” position and the centrifuge should be restarted by pressing the start (green) button.
9. An appropriate UPS may be used for un - interrupted function.

Specifications:

Choice of two speeds (low: 6,000 rpm and high: 8,000 rpm).

Adjustable timer in steps of 15 seconds (15 seconds to infinity).

Safety switch turns the centrifuge "off " on opening the lid.